



PRODUCT BULLETIN


SEPOR, INC.
718 N. FRIES AVE.
P. O. BOX 578
WILMINGTON, CA 90748
PHONE: 310-830-6601
FAX: 310-830-9336

Web Site: <http://www.sepor.com>

Automagnet Operation and Maintenance

Pick up automagnet, placing fingers under top of body, place magnet in material and depress plunger with thumb. Holding plunger down, move bottom of magnet around in material. Magnetic particles will stick to outside of automagnet body, on bottom of body. Remove automagnet from material and move to clean container, slowly release plunger and magnet material will fall from plunger. Repeat, as required.

In the event the plunger does not depress, or has difficulty returning to the release position, follow instructions below for repair of abused units.

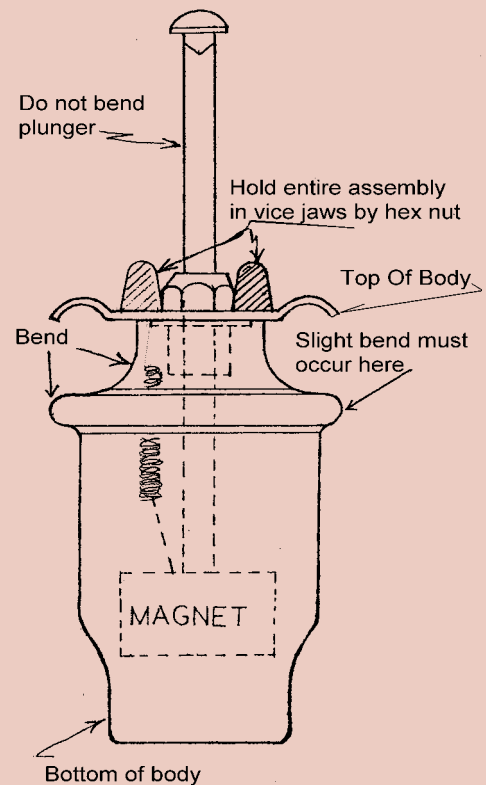

SEPOR AUTOMAGNET

REPAIR OF ABUSED OR DROPPED UNITS

1. Lift Automagnet to eye level and sight the plunger angle-incline.
2. Prepare to bend top of body in opposite direction of plunger-incline.
3. Insert Automagnet in vice jaws and hold by the hex nut. **DO NOT PUT VICE JAWS ON THE PLUNGER.**
4. Grasp the bottom of the body with the palm of the hand and bend top of body (slightly and gently) to make plunger in line with the main body, or so that the interior magnet assembly does not touch the sides of the bottom section.
5. SEPOR, Inc. will not accept responsibility for any adjustments made by the operator.